### Practical Session N°4: Observation of Cellular Organelles under the Microscope

# **Objectives**

- Learn how to prepare slides for microscopic observation.
- Identify the main cellular organelles visible under an optical microscope.
- Compare the cellular structures of plant and animal cells.

# **Required Materials**

- 1. Optical microscope.
- 2. Slides and coverslips.
- 3. Biological stains (methylene blue, methyl green, iodine solution, neutral red).
- 4. Pasteur pipettes.
- 5. Distilled water.
- 6. Filter paper or blotting paper.
- 7. Scalpel or razor blade.
- 8. Fine forceps.
- 9. Specimens for observation:
  - o **Plant cells**: onion skin, epidermis of Elodea leaf or other aquatic plants.
  - Animal cells: buccal cell smear.
  - Yeast cells.

#### **Procedure**

### 1. Observation of the Nucleus (Animal Cells)

## • Preparation:

- o Gently collect buccal cells using a cotton swab or a clean slide edge.
- o Place the cells on a slide and add one drop of methylene blue.
- o Carefully cover with a coverslip, avoiding air bubbles.

#### Observation:

- o Observe under the microscope at low magnification (10x), then increase to 40x.
- o Identify the nucleus, which will appear dark blue.

### 2. Observation of Chloroplasts (Plant Cells)

# • Preparation:

- o Take a small piece of Elodea leaf or onion skin.
- o Place the fragment on a slide with a drop of distilled water.

o Cover with a coverslip and blot excess water with filter paper.

#### • Observation:

o Observe under the microscope. Chloroplasts appear as green oval structures.

## 3. Observation of Mitochondria (Special Staining)

### • Preparation:

- o Use yeast cells or buccal cells.
- o Add a drop of neutral red, a stain that highlights mitochondria.
- o Cover with a coverslip.

## • Observation:

o Look for small red dots dispersed in the cytoplasm.

### 4. Observation of Vacuoles (Plant Cells)

# • Preparation:

- Prepare a slide with a piece of onion skin and add a drop of iodine solution (Lugol's iodine).
- o Cover with a coverslip.

### • Observation:

 Vacuoles appear as clear structures in the cytoplasm, surrounded by stained cell walls and nuclei.

# **Analysis and Interpretation**

- 1. Draw annotated diagrams for each type of cell observed.
- 2. Compare the structures of animal and plant cells (e.g., presence of a cell wall, chloroplasts, and vacuole in plant cells; absence of these organelles in animal cells).